

P30XL Tip System Technique Modification for Biomek FX and NX systems

After installing the P30XL tip definition, the existing system techniques will require modification to allow use of the new tip.

For Biomek FX systems:

1. Open the Technique Browser
2. Select *1536 Plate*
3. Click on Properties
4. Scroll to the tip section at the bottom of the list
5. Select P30XL
6. Change the Maximum Volume to 50
7. Select O.K.
 - Select the *Low Volume 384 Plate* and repeat steps 3 to 7.
 - Select the *Washing* technique and repeat steps 3 to 5.

For Biomek NX systems:

1. Open the Technique Browser
2. Select *1536 Plate*
3. Click on Properties
4. Scroll to the tip section at the bottom of the list
5. Select P30XL
6. Change the Maximum Volume to 50
7. Select O.K.
 - Select *Default-AP384* and repeat steps 3 to 5.
 - Select *Washing 384* and repeat steps 3 to 5.
 - Select *AP384-Low Volume* and repeat steps 3 to 5.